

# GENLISA™ PLANT DIAGNOSTIC DAS ELISA

**REF** : DASREAG

Ver 1.0

**RUO**

<b>RUO</b>	For Research & Use Only	<b>REF</b>	Catalog Number
	Store At	<b>LOT</b>	Batch Code
	Manufactured By		Biological Risk
	Expiry Date		Consult Operating Instructions

*For Research Use Only. Purchase does not include or carry the right to resell or transfer this product either as a stand-alone product or as a component of another product. Any use of this product other than the permitted use without the express written authorization of KRISHGEN BioSystems is strictly prohibited.*



**KRISHGEN BioSystems**

Unit Nos#318/319, Shah & Nahar,  
Off Dr E Moses Road, Worli, Mumbai 400018.  
Tel: 91 (22) 49198700 | Email: sales@krishgen.com

## ELISA Reagents DAS ELISA, alkaline phosphatase conjugate

### Components Available in the Kit:

Pack Size	500 assays	1000 assays	5000 assays	Storage
Coating antibody	0.25 ml	0.50 ml	2 x 1.25 ml	4°C
Detecting conjugate alkaline phosphatase	0.25 ml	0.50 ml	2 x 1.25 ml	4°C
96 well ELISA plates	5	10	50	Room Temperature
Instruction Manual	1	1	1	

### Safety and Storage:

Always wash hands thoroughly after using this product. Prevent direct skin and eye contact with, or ingestion of product components. Obtain medical attention in case of accidental ingestion of reagent components.

All reagent components should be stored at the recommended temperature to assure their full shelf life. Do not store prepared working solution from day to day.

Please contact Krishgen Biosystems, if you have any questions about safety and storage of this product.

### Preparing For The Test:

- 1) Check all the components in the package of ELISA Reagents.
- 2) Prepare all buffer solutions according to the attached buffer formulations.
- 3) Make sure all laboratory equipment and facilities required for the test are ready.
- 4) Prepare a humid box for incubation steps.
- 5) Make a copy of the attached recording sheet and create a loading diagram by recording the locations of your samples, controls, and other reagents needed.

### Coating Plate with Antibody:

Lay out all items that will be required for the plate coating step before beginning. Prepare coating antibody in a container made of glass, polyethylene or any material that does not readily bind coating antibody. Coat the plate immediately after preparing the coating antibody. Some coating antibody can be lost if too much time elapses between diluting the coating

### Antibody and Coating the Plate:

- 1) The volume of coating buffer required depends on the number of test wells used; 100 ul is needed per test well. One way to estimate the volume needed is to prepare 1 ml of coating buffer for each 8 well strip used, or 10 ml for each 96 well plate.
- 2) Dilute the concentrated coating antibody into coating buffer at the dilution given on the label. Mix well. Always prepare coating antibody immediately before use.

- 3) Pipette 100 ul of coating antibody into each well.
- 4) Incubate the plate in a humid box for overnight in the refrigerator (4°C) or 4 hours at room temperature (21-24°C).

**Preparing Samples:**

- 1) Select symptomatic and/or infective tissues for the test. Leaf tissue is often used in ELISA testing. Plant tissues such as stem, sprout, seed, tuber, root and others can also be used.
- 2) We suggest that each test well be used for only one sample. In some cases, composites of up to ten leaves per test well can be used to make testing more economical. However, too many plant samples per well can reduce the sensitivity of the test.
- 3) KRISHGEN's KBBR01 buffer can be used as extraction buffer for most of the plant samples. However, other buffers are also recommended for some plant species.
- 4) Grind sample with a mortar and pestle, or other grinding device. If you are using a mortar and pestle, wash and rinse it thoroughly between samples.
- 5) If you extract plant sap, dilute the sap into sample extraction buffer at a ratio of 1:10 (sap volume : buffer volume) or you can grind plant tissue in extraction buffer at a 1:10 ratio (tissue weight : Extraction Buffer volume).
- 6) If you have any questions about sampling, sample preparation, or the appropriate extraction buffer for your samples, please contact Krishgen Biosystems.

**Plate Washing:**

- 1) Wash the plate when the incubation is complete. Use a quick flipping motion to empty the wells into a sink or waste container.
- 2) Wash the plate by filling the wells with PBST, then quickly emptying them again. Repeat 4 to 6 times.
- 3) To remove drops of PBST from the wells after washing, hold the frame upside down and tap firmly on a folded paper towel.

**Sample Dispensing and Incubation:**

About 100 ul of diluted sample extract is needed per test well. Always have an additional amount to assure easy dispensing. A convenient way to prepare this diluted sample is to measure 100 ul of undiluted sap into a small test tube and then add 1 ml of extraction buffer.

Following your loading diagram on your recording sheet, dispense 100 ul of prepared sample into sample wells. Dispense 100 ul of positive control into positive control wells, and dispense 100 ul of negative control or extraction buffer into negative control wells.

Put the plate inside the humid box and incubate for 2.5 hours at room temperature (21-24°C) or overnight in the refrigerator (4°C).

**Preparing Enzyme Conjugate:**

Always make enzyme conjugate solution within 10 minutes before use. Prepare the enzyme conjugate, using KRISHGEN's KBCONJ01 buffer and a cleaning container.

The volume of KBCONJ01 buffer required depends on the number of test wells used; 100 ul are needed per test well. To estimate the volume needed, prepare 1 ml for each 8 well strip used, or 10 ml for each 96 well plate.

The volume of enzyme conjugate required for each test is calculated based on the volume of KBCONJ01 buffer used and on the dilutions given on the bottles. Use a new, sterile pipette tip and change the tip for each pipetting to prevent contamination.

First dispense appropriate volume of KBCONJ01 buffer into a cleaning container and then add enzyme conjugate according to the dilution given on the label. Mix the conjugate solution thoroughly. It is important to mix the enzyme conjugate well for a consistent test result.

Prepare enzyme conjugate just before use. Keep the prepared enzyme conjugate at a safe place and use it after washing the plate.

**Washing Plate:**

Wash the plate when the incubation is complete. Use a quick flipping motion to empty the wells into a sink or waste container without mixing the contents.

Wash the plate by filling the wells with PBST, then quickly emptying them again. Repeat 6 to 8 times. To remove drops of PBST from the wells after washing, hold the frame upside down and tap firmly on a folded paper towel.

**Enzyme Conjugate Incubation:**

Dispense 100 ul of prepared enzyme conjugate per well for all test wells. Incubate the plate in the humid box for 2.5 hours at room temperature (21-24°C).

**Preparing Substrate Solution:**

Concentration of PNP in substrate is 1 mg/ml. Each PNP tablet will make 5 ml of PNP solution, which is enough for 48 test wells or five 8 well strips.

Do not touch the PNP tablets or expose the PNP solution to strong light. Light or contamination could cause background color in negative wells.

Prepare PNP substrate about 10-15 minutes before the end of the above incubation step. Measure 5 ml of PNP buffer for each tablet and then add the PNP tablets to the buffer. Mix by vortexing or stirring to let the PNP tablet fully dissolve in the buffer.

**Washing the Plate:**

Wash the plate 6 to 8 times with PBST as instructed above.

**Incubation with Substrate:**

Dispense 100 ul of PNP substrate solution per well. Incubate the plate for 30 to 60 minutes in a humid box at room temperature (21-24°C).

To stop reaction, add 50 ul of 3M Sodium Hydroxide to each well. This step is optional. The plate can be interpreted visually or with a plate reader without adding the Stop Solution.

**Data Interpretation:**

Test results can be examined by eye, or measured on a plate reader at 405 nm.

Development of yellow color in test wells indicates positive results. Wells in which there is no significant color development indicate negative results. Test results are valid only if positive control wells give a positive result and negative control wells remain clear.

Results may be interpreted after more than 60 minutes of incubation as long as negative control wells remain virtually clear.

**Buffer Formulations****Coating Buffer**

Sodium Carbonate (anhydrous)	1.60 gm
Sodium Bicarbonate	2.92 gm
Sodium Azide	0.20 gm

Dissolve in distilled water and make to 1000 ml. Adjust pH to 9.6. Store at 4°C.

**PBST Buffer**

Sodium Phosphate, dibasic, (anhydrous)	1.15 gm
Potassium Phosphate, monobasic (anhydrous)	0.20 gm
Sodium Chloride	8.00 gm
Potassium Chloride	0.20 gm
Tween-20	0.50 gm

Dissolve in distilled water and make to 1000 ml. Adjust pH to 7.3.

**KBBR01 Buffer**

Powdered Egg (chicken) albumin, Grade II	2.00 gm
Polyvinyl Pyrrolidone (PVP) MW24-40,000	10.00 gm
Sodium Sulfite (anhydrous)	1.30 gm
Sodium Azide	0.20 gm
Tween-20	10.00 gm

Dissolve in 1000 ml of 1X PBST. Adjust pH to 7.3. Store at 4°C.

**KBCONJ01 Buffer**

Bovine Serum Albumin (BSA)	2.00 gm
Polyvinyl Pyrrolidone (PVP) MW24-40,000	10.00 gm
Sodium Azide	0.20 gm

Dissolve in 1000 ml of 1X PBST. Adjust pH to 7.3. Store at 4°C.

**PNP Buffer**

Diethanolamine	97.00 ml
Magnesium Chloride	0.10 gm
Sodium Azide	0.20 gm

Dissolve in 800 ml distilled water. Adjust pH to 9.8 with hydrochloric acid. Adjust final volume to 1000 ml with distilled water. Store at 4° C.

**RECORDING SHEET FOR ELISA**

TEST: \_\_\_\_\_  
DATE: \_\_\_\_\_  
BY: \_\_\_\_\_

TIMING:  
Coating - \_\_\_\_\_  
Sample - \_\_\_\_\_  
EC - \_\_\_\_\_  
Substrate - \_\_\_\_\_

KEY POINTS:  
Coating Antibody: \_\_\_\_\_ ul  
Coating Buffer: \_\_\_\_\_ ml  
Enzyme Conjugate: \_\_\_\_\_ ul  
KBCONJ1: \_\_\_\_\_ ml  
PNP Substrate: \_\_\_\_\_ ml

1      2      3      4      5      6      7      8      9      10      11      12

A												
B												
C												
D												
E												
F												
G												
H												

**RESULTS/CONCLUSIONS:**

- 1. \_\_\_\_\_
- 2. \_\_\_\_\_
- 3. \_\_\_\_\_

**LIMITED WARRANTY**

Krishgen Biosystems does not warrant against damages or defects arising in shipping or handling, or out of accident or improper or abnormal use of the Products; against defects in products or components not manufactured by Krishgen Biosystems, or against damages resulting from such non-Krishgen Biosystems made products or components. Krishgen Biosystems passes on to customer the warranty it received (if any)

from the maker thereof of such non Krishgen made products or components. This warranty also does not apply to Products to which changes or modifications have been made or attempted by persons other than pursuant to written authorization by Krishgen Biosystems.

THIS WARRANTY IS EXCLUSIVE. The sole and exclusive obligation of Krishgen Biosystems shall be to repair or replace the defective Products in the manner and for the period provided above. Krishgen Biosystems shall not have any other obligation with respect to the Products or any part thereof, whether based on contract, tort, and strict liability or otherwise. Under no circumstances, whether based on this Limited Warranty or otherwise, shall Krishgen Biosystems be liable for incidental, special, or consequential damages.

This Limited Warranty states the entire obligation of Krishgen Biosystems with respect to the Products. If any part of this Limited Warranty is determined to be void or illegal, the remainder shall remain in full force and effect.

**Krishgen Biosystems. 2018**

**THANK YOU FOR USING KRISHGEN PRODUCT!**

*Note: GENLISA™ is a registered trademark of KRISHGEN BIOSYSTEMS.*