






KRISHZYME™ Swift™ PNGase-F, lyophilized

REF : KPGF-001

Ver 1.0

RIUO

RIUO	For Research & Industrial Use Only	REF	Catalog Number
	Store At	LOT	Batch Code
	Manufactured By		Biological Risk
	Expiry Date		Consult Operating Instructions

For Research and Industrial Use Only. Purchase does not include or carry the right to resell or transfer this product either as a stand-alone product or as a component of another product. Any use of this product other than the permitted use without the express written authorization of KRISHGEN BioSystems is strictly prohibited.

REF KPGF-001

KRISHGEN BioSystems For US/Europe Customers: toll free +1(888)-970-0827 | tel +1(562)-568-5005
For Asia/India Customers: +91(22)-49198700
Email: sales1@krishgen.com | <http://www.krishgen.biz>

Peptide N-glycosidase F, commonly referred to as PNGase F, is an amidase of the peptide-N4-(N-acetyl-beta-glucosaminyl) asparagine amidase class. PNGase F is the most effective enzymatic method for removing almost all N-linked oligosaccharides from glycoproteins.

High Mannose oligosaccharide

Hybrid oligosaccharide

Complex oligosaccharides

Asn

GlcNAc

Mannose

Galactose

Sialic

Fucose

Product Size :

Catalog number	Pack Size	Concentration
KPGF-001-A	50,000U / 50 ul	1,000,000 U /ml
KPGF-001-B	2 x 50,000U / 50 ul	

KRISHZYME™ Swift™ PNGase F is supplied as a vial of enzyme containing 100 ug of PNGase-F lyophilized from 20mM Tris-HCl, 50mM NaCl, pH 7.5. Resuspend in 50 ul of double distilled water to get a concentration of 1,000 U/ul or 50,000 U/vial.

The following reagents are supplied with this product:

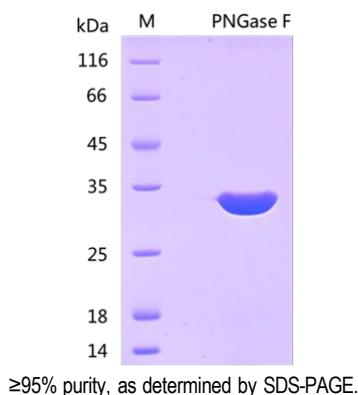
Composition	Formula	Concentration
Denaturing Buffer	5%SDS, 0.4M DTT	10X
Assay Buffer 2	0.5M Sodium Phosphate (pH7.5 at 25°C)	10X
NP-40 Solution		10%

Product Source:

Recombinant gene cloned from *Flavobacterium meningsepticum* and expressed in *E. coli*.

Product Quality:

≥95% purity, as determined by SDS-PAGE. No other exoglycosidase, endoglycosidase, and protease activity were contaminated.

**Unit Definition:**

One unit of PNGase F is defined as the amount of enzyme required to remove > 95% of the carbohydrate from 10 ug of denatured RNase B in 1 hour at 37°C in a total reaction volume of 10 ul.

Storage Temperature:

4°C

Characteristic:

- Recombinant enzyme
- Glycerol-free for optimal performance in HPLC and mass spectrometry analysis
- ≥95% purity, as determined by SDS-PAGE
- Optimal activity and stability for up to 12 months
- Can be used under native or denaturing conditions
- Optimized for deglycosylation of glycoproteins; leaves N-glycan core oligosaccharides intact and suitable for further analysis

Applications:

- Characterizing whether the protein is glycosylated
- Determining the location of glycosylation on the protein
- Characterizing the glycan structure
- Protein trafficking
- Release of intact N-linked glycans from glycopeptides and glycoproteins
- Structure-function studies of N-glycosylated glycoproteins
- Preparation of deglycosylated proteins for molecular weight estimation or crystallography studies

Suggestions for Use:**Denaturing Reaction Conditions:**

- 1) Combine 10 - 100 ug of glycoprotein, 1 ul of 10×Denaturing Buffer and H₂O (if necessary) to make a 10 ul total reaction volume;
- 2) Denature glycoprotein by heating reaction at 100°C for 10 minutes;
- 3) Chill denatured glycoprotein on ice and centrifuge 10 seconds;
- 4) Make a total reaction volume of 20 ul by adding 2 ul 10×Assay Buffer 2, 2 ul 10% NP- 40 Solution and 6 ul H₂O;
- 5) Add 1 ul PNGase F, mix gently;
- 6) Incubate reaction at 37°C for 1-3 hours.
- 7) Analyze by method of choice.

Non-Denaturing Reaction Conditions:

- 1) Combine 10 - 100 ug of glycoprotein, 2 ul of (10X) Assay Buffer 2, 2-5 ul PNGase F (Glycerol-free) and H₂O (if necessary) to make a 20 ul total reaction volume, mix gently;
- 2) Incubate reaction at 37°C for 4-24 hours;
- 3) Analyze by method of choice.

Notes :

- When deglycosylating a native glycoprotein it is recommended that an aliquot of the glycoprotein is subjected to the denaturing protocol to provide a positive control for the fully deglycosylated protein. The non-denatured reaction can then be compared to the denatured reaction to determine the extent of reaction completion;
- To deglycosylate a native glycoprotein, longer incubation time as well as more enzyme may be required;
- The simplest method of assessing the extent of deglycosylation is by mobility shifts on SDS-PAGE gels;
- Since PNGase F (Glycerol-free), Recombinant activity is inhibited by SDS, it is essential to have NP-40 in the reaction mixture. It is not known why this non-ionic detergent counteracts the SDS inhibition at the present time;
- PNGase F (Glycerol-free), Recombinant will not cleave N-linked glycans containing core α1-3 Fucose;
- Recommended storage temperature is 4°C, avoid repeat freeze-thaw cycles.

References:

Characterization of glycoproteins and their associated oligosaccharides through the use of endoglycosidases. ... F Maley, RB Trimble, AL Tarentino... - Analytical biochemistry, 1989 - Elsevier

Deglycosylation of asparagine-linked glycans by peptide: N-glycosidase F. ... AL Tarentino, CM Gomez, TH Plummer Jr - Biochemistry, 1985 - ACS Publications/

Demonstration of peptide: N-glycosidase F activity in endo-beta-N-acetylglucosaminidase F preparations. ... TH Plummer, JH Elder, S Alexander, AW Phelan... - Journal of Biological ..., 1984 - ASBMB

Glycosylation of Pichia pastoris -derived proteins. ... RK Bretthauer, FJ Castellino - Biotechnology and applied ..., 1999 - Wiley Online Library.

Pharmacological chaperones rescue cell-surface expression and function of misfolded V2 vasopressin receptor mutants. ... JP Morello, A Salahpour, A Laperrière... - The Journal of ..., 2000 - Am Soc Clin Investig.

A monolithic PNGase F enzyme microreactor enabling glycan mass mapping of glycoproteins by mass spectrometry. ... AK Palm, MV Novotny - ... An International Journal Devoted to the ..., 2005 - Wiley Online Library.

Protein-protein interaction and not glycosylation determines the binding selectivity of heterodimers between the calcitonin receptor-like receptor and the ... S Hilaiet, SM Foord, FH Marshall, M Bouvier - Journal of Biological ..., 2001 - ASBMB

Purification and Structure-Function Analysis of Native, PNGase F-Treated, and Endo-. beta.-galactosidase-Treated CHIP28 Water Channels. AN Van Hoek, MC Wiener, JM Verbavatz, D Brown... - Biochemistry, 1995 - ACS Publications

Isolation and characterization of CD47 glycoprotein: a multispinning membrane protein which is the same as integrin-associated protein (IAP) and the ovarian tumour ... WJ Mawby, CH Holmes, DJ Anstee, FA Spring... - Biochemical ..., 1994 - biochemj.org

DC-SIGN and L-SIGN are high affinity binding receptors for hepatitis C virus glycoprotein E2 ..., S Fong, A Amara, C Houles, F Fieschi... - Journal of Biological ..., 2003 - ASBMB

LIMITED WARRANTY

Krishgen Biosystems does not warrant against damages or defects arising in shipping or handling, or out of accident or improper or abnormal use of the Products; against defects in products or components not manufactured by Krishgen Biosystems, or against damages resulting from such non-Krishgen Biosystems made products or components. Krishgen Biosystems passes on to customer the warranty it received (if any) from the maker thereof of such non Krishgen made products or components. This warranty also does not apply to Products to which changes or modifications have been made or attempted by persons other than pursuant to written authorization by Krishgen Biosystems.

THIS WARRANTY IS EXCLUSIVE. The sole and exclusive obligation of Krishgen Biosystems shall be to repair or replace the defective Products in the manner and for the period provided above. Krishgen Biosystems shall not have any other obligation with respect to the Products or any part thereof, whether based on contract, tort, and strict liability or otherwise. Under no circumstances, whether based on this Limited Warranty or otherwise, shall Krishgen Biosystems be liable for incidental, special, or consequential damages.

This Limited Warranty states the entire obligation of Krishgen Biosystems with respect to the Products. If any part of this Limited Warranty is determined to be void or illegal, the remainder shall remain in full force and effect.

Krishgen Biosystems. 2025

THANK YOU FOR USING KRISHGEN PRODUCT!

KRISHGEN BIOSYSTEMS®, GENLISA®, DHARMAPLEX™, GENBULK™, GENLISA™, KRISHZYME®, KRISHZYME SIWFT®, KRISHGEN®, KRIBIOLISA®, KRISHPLEX®, TITANIUM®, QUALICHEK® are registered trademarks of KRISHGEN BIOSYSTEMS PRIVATE LIMITED. ©KRISHGEN BIOSYSTEMS. ALL RIGHTS RESERVED.

KRISHGEN BIOSYSTEMS | OUR REAGENTS | YOUR RESEARCH |