






# KRISHZYME™ $\alpha$ 2-3,6,8,9 Neuraminidase

**REF** : KPGF-005

Ver 2.0

**RIUO**

<b>RIUO</b>	<b>For Research &amp; Industrial Use Only</b>	<b>REF</b>	<b>Catalog Number</b>
	<b>Store At</b>	<b>LOT</b>	<b>Batch Code</b>
	<b>Manufactured By</b>		<b>Biological Risk</b>
	<b>Expiry Date</b>		<b>Consult Operating Instructions</b>

*For Research and Industrial Use Only. Purchase does not include or carry the right to resell or transfer this product either as a stand-alone product or as a component of another product. Any use of this product other than the permitted use without the express written authorization of KRISHGEN BioSystems is strictly prohibited.*


**REF** KPGF-004

**KRISHGEN BioSystems** For US/Europe Customers: toll free +1(888)-970-0827 | tel +1(562)-568-5005  
For Asia/India Customers: +91(22)-49198700  
Email: sales1@krishgen.com | <http://www.krishgen.biz>

### Product Description:

Neuraminidase is the common name for Acetyl-neuraminyl hydrolase (Sialidase).  $\alpha$ 2-3,6,8,9 Neuraminidase is a broad specificity sialidase, which cleaves linear and branched non-reducing terminal sialic acid residues from glycoproteins, glycopeptides, and oligosaccharides. It can be used for glycan analysis and characterization and intact glycoprotein remodeling, in vitro and in vivo.

KRISHZYME™  $\alpha$ 2-3,6,8,9 Neuraminidase is a recombinant glycosidase cloned from *Arthrobacter ureafaciens* and expressed in *E. coli*. The enzyme releases  $\alpha$ (2,3)-,  $\alpha$ (2,6)-,  $\alpha$ (2,8)-, and  $\alpha$ (2,9)-linked N-acetylneuraminic acid from complex carbohydrates. The initial rate of hydrolysis of  $\alpha$ (2,6) linkages is reported to be approximately twice that of  $\alpha$ (2,3)-linked sialic acid, however, in practice, this kinetic selectivity is of little consequence during extended incubations. Effective digestion of glycolipid substrates is facilitated by addition of a detergent, such as sodium taurodeoxycholate to the incubation.

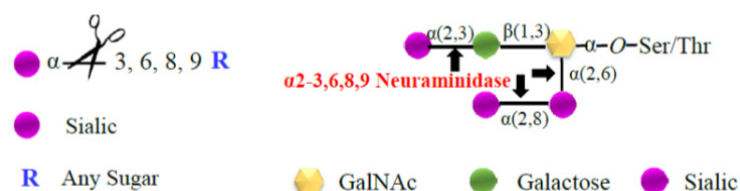

The enzyme releases  $\alpha$ (2,3)-,  $\alpha$ (2,6)-,  $\alpha$ (2,8)-, and  $\alpha$ (2,9)-linked N-acetylneuraminic acid from complex carbohydrates.

### Product Size:

Catalog Number	Pack Size	Concentration
KPGF-005-A	0.6 U / 30 ul	20 U /ml
KPGF-005-B	5 x 0.6 U / 30 ul	

### Physical Form:

KRISHZYME™  $\alpha$ 2-3,6,8,9 Neuraminidase is supplied as a liquid in 20mM Tris-HCl (pH 7.5 at 25°C), 50 mM NaCl, 1mM EDTA at a concentration of 20 U/ml.

### Reagents Supplied:

The following reagents are supplied with this product:

Composition	Formula	Concentration
Assay Buffer 1	50 mM CaCl <sub>2</sub> , 500 mM sodium acetate, pH 5.5 at 25°C	10X

### Product Source:

Recombinant gene cloned from *Arthrobacter ureafaciens* and expressed in *E. coli*. Krishzyme™  $\alpha$ 2-3,6,8,9 Neuraminidase has a molecular weight of 66kDa.

### Product Quality:

≥95% purity, as determined by SDS-PAGE. No other exoglycosidase, endoglycosidase, and protease activity were contaminated.

**Unit Definition:**

One unit is defined as the amount of enzyme required to catalyze the release of 1 umole of p-nitrophenol from p-nitrophenyl- $\alpha$ -D-N-acetylneuraminic acid per minute at 37° C, pH 5.5.

**Storage Temperature:**

Store at 4°C, in a frost free refrigerator

**Characteristic:**

- Recombinant enzyme with no detectable endoglycosidase or other exoglycosidases contaminating activities
- $\geq 95\%$  purity, as determined by SDS-PAGE
- Optimal activity and stability for up to 24 months
- Can be used under native or denaturing conditions
- Acts on both Neu5Ac and Neu5Gc
- Double digest with other exoglycosidases and endoglycosidases
- Tolerant of moderate levels (0.5-1.0%) of detergents

**Applications:**

- Structural analysis of oligosaccharides
- Determining sialic acid linkage
- Glycoprotein deglycosylation
- Removing heterogeneity from glycoproteins

**Suggestions for Use:**

- 1) Combine 1-100 ug of glycoprotein and H<sub>2</sub>O (if necessary) in a total reaction volume of 8 ul;
- 2) Add 1 ul of 10X Assay Buffer 1 to make a 10 ul total reaction volume;
- 3) Add 1 ul alpha2-3,6,8,9 Neuraminidase, mix gently ;
- 4) Incubate at 37°C for 1 hour.

**Notes :**

- The amount of exoglycosidase enzyme required varies when different substrates are used. Start with 1-2 ul for 1-100ug of glycoprotein for one hour in a 10-25  $\mu$ l reaction. If there is still undigested material, let the reaction go overnight;
- The reaction can be scaled up linearly;
- Higher concentrations of enzyme as well as longer incubation times may be necessary for cleavage of branched structures.

**References:**

Characterization of glycoproteins and their associated oligosaccharides through the use of endoglycosidases. ... F Maley, RB Trimble, AL Tarentino... - Analytical biochemistry, 1989 - Elsevier

Deglycosylation of asparagine-linked glycans by peptide: N-glycosidase F. ... AL Tarentino, CM Gomez, TH Plummer Jr - Biochemistry, 1985 - ACS Publications/

Demonstration of peptide: N-glycosidase F activity in endo-beta-N-acetylglucosaminidase F preparations. ... TH Plummer, JH Elder, S Alexander, AW Phelan... - Journal of Biological ..., 1984 - ASBMB

Glycosylation of *Pichia pastoris* -derived proteins. ... RK Bretthauer, FJ Castellino - Biotechnology and applied ..., 1999 - Wiley Online Library.

Pharmacological chaperones rescue cell-surface expression and function of misfolded V2 vasopressin receptor mutants. ... JP Morello, A Salahpour, A Laperrière... - The Journal of ..., 2000 - Am Soc Clin Investig.

A monolithic PNGase F enzyme microreactor enabling glycan mass mapping of glycoproteins by mass spectrometry. ... AK Palm, MV Novotny - ... An International Journal Devoted to the ..., 2005 - Wiley Online Library.

Protein-protein interaction and not glycosylation determines the binding selectivity of heterodimers between the calcitonin receptor-like receptor and the ... S Hilairiet, SM Foord, FH Marshall, M Bouvier - Journal of Biological ..., 2001 - ASBMB

Purification and Structure-Function Analysis of Native, PNGase F-Treated, and Endo-. beta.-galactosidase-Treated CHIP28 Water Channels. AN Van Hoek, MC Wiener, JM Verbavatz, D Brown... - Biochemistry, 1995 - ACS Publications

Isolation and characterization of CD47 glycoprotein: a multispanning membrane protein which is the same as integrin-associated protein (IAP) and the ovarian tumour ... WJ Mawby, CH Holmes, DJ Anstee, FA Spring... - Biochemical ..., 1994 - biochemj.org

## LIMITED WARRANTY

Krishgen Biosystems does not warrant against damages or defects arising in shipping or handling, or out of accident or improper or abnormal use of the Products; against defects in products or components not manufactured by Krishgen Biosystems, or against damages resulting from such non-Krishgen Biosystems made products or components. Krishgen Biosystems passes on to customer the warranty it received (if any) from the maker thereof of such non Krishgen made products or components. This warranty also does not apply to Products to which changes or modifications have been made or attempted by persons other than pursuant to written authorization by Krishgen Biosystems.

THIS WARRANTY IS EXCLUSIVE. The sole and exclusive obligation of Krishgen Biosystems shall be to repair or replace the defective Products in the manner and for the period provided above. Krishgen Biosystems shall not have any other obligation with respect to the Products or any part thereof, whether based on contract, tort, and strict liability or otherwise. Under no circumstances, whether based on this Limited Warranty or otherwise, shall Krishgen Biosystems be liable for incidental, special, or consequential damages.

This Limited Warranty states the entire obligation of Krishgen Biosystems with respect to the Products. If any part of this Limited Warranty is determined to be void or illegal, the remainder shall remain in full force and effect.

**Krishgen Biosystems. 2025**

## THANK YOU FOR USING KRISHGEN PRODUCT!

KRISHGEN BIOSYSTEMS®, GENLISA®, DHARMAPLEX™, GENBULK™, GENLISA™, KRISHZYME®, KRISHZYME SIWFT®, KRISHGEN®, KRIBIOLISA®, KRISHPLEX®, TITANIUM®, QUALICHEK® are registered trademarks of KRISHGEN BIOSYSTEMS PRIVATE LIMITED. ©KRISHGEN BIOSYSTEMS. ALL RIGHTS RESERVED.

**KRISHGEN BIOSYSTEMS | OUR REAGENTS | YOUR RESEARCH |**